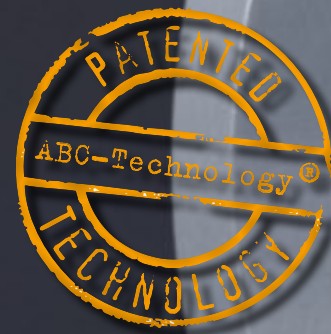
## Anti-Bearding-Technik<sup>®</sup> speziell für Pharma & Food *Anti-Bearding Technology<sup>®</sup> especially for Pharma & Food*

- Kein Produktaufbau dank SCHLICK Anti-Bearding-Cap (ABC)  
*No product build-up thanks to the SCHLICK Anti-Bearding-Cap (ABC)*

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# Anti-Bearding for Pharma and Food 100% made in Germany

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*Optimal homogenous and reproducible spray results*

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### SCHLICK nozzle technology increases productivity

The newly developed SCHLICK Anti-Bearding Technology® offers optimum operating security within coating-facilities. The patented air cap for flat spray two-substance nozzles, the so-called „Anti-Bearding Cap“ (ABC), crucially prevents turbulence in the area of the air cap and the nozzle exit, by means of its novel rounded shape.

### Perfect 'beardless' spraying

Build-up, caking, „bearding“ or clogging of the nozzle exit are prevented as far as possible in this way. Interruptions to the coating process, as a consequence of cleaning work on the nozzle, are almost completely omitted. The SCHLICK ABC spray features an exceedingly homogenous liquid distribution, in a simultaneously very fine, uniform and reproducible drop size distribution.

The SCHLICK Anti-Bearding Technology® is provided for lab coaters and for production facilities. A simple scale up is possible.

Modell	Nano-Düse 951 7-1 S24 NANO ABC	Labordüse 970 7-1 S75 ABC	Produktionsdüse 930 7-1 S35 ABC	Produktionsdüse 930 7-1 S45 ABC
Model	Nano nozzle 951 7-1 S24 NANO ABC	Laboratory nozzle 970 7-1 S75 ABC	Production nozzle 930 7-1 S35 ABC	Production nozzle 930 7-1 S45 ABC
Trommelgröße Drum size	8,5" / 10,5" / 12"	12" / 15" / 19"	19" / 24" und größer 19" / 24" and above	19" / 24" und größer 19" / 24" and above
Streukegel Spray angle	10 – 60°	10 – 60°	10 – 60°	60°
Zerstäubungsform Spray pattern	Ellipsenförmiger Flachstrahl Oval flat spray	Ellipsenförmiger Flachstrahl Oval flat spray	Ellipsenförmiger Flachstrahl Oval flat spray	Ellipsenförmiger Flachstrahl Oval flat spray
Durchsatz Capacity	1 – 10 g/min (Coating)	5 – 60 g/min (Coating)	30 – 180 g/min (Coating)	30 – 180 g/min (Coating)

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**Innovative product design**

- GMP design: Simple construction (7 individual parts + 4 O rings)
- Cleaning needle
- Liquid return system is available
- Manual assembly / disassembly
- Easy cleaning
- Materials conform to FDA: 1.4404 (AISI 316 L), EPDM-O-rings

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Flow behaviour with conventional flat spray cap and the consequence to the bearding effect

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Flow behaviour with SCHLICK Anti-Bearding Cap (ABC) and the consequence to the bearding effect

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A comparison of the flow patterns reveals that the dust particle flow (simulated here in the form of fog) stays away from the anti-bearding cap better than it does from the conventional flat spray cap.

- Produktübersicht  
Product Overview
- Zweistoffdüsen  
Two-Substance  
Nozzles
- ABC/PCA-1 Technik  
ABC/PCA Technique
- Drei-/Vier-Stoffdüsen  
Three-/Four-  
Substance Nozzles
- Multispray  
Multispray
- Einsteckrohre  
Insertion Pipes
- Düsenköpfe  
Nozzle Heads
- Vollkegeldüsen  
Full Cone Nozzles
- Hohlkegeldüsen  
Hollow Cone Nozzles
- Hartmetaldüsen  
Carbide Nozzles
- Flachstrahldüsen  
Flat Spray Nozzles
- Glattdüsen  
Smooth-Jet Nozzles
- Mischdüsen  
Mixing Nozzles
- Dampf-/  
Luftstrahldüsen  
Laval Nozzles
- Reinigungsdüsen  
Cleaning Nozzles
- Zubehör  
Accessories
- Test Center  
Test Center

# Anti-Bearding-Technik® im Detail

## Anti-Bearding Technology® in Detail

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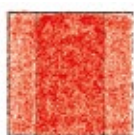
### ABC spray: perfect droplets in the laboratory and production

The ideal ABC spray is characterised by a very even ellipsis. This is achieved by means of a balance between the pressures – the forces of the **atomising air (AA)** and the **pattern air (PA)**.

In the following figure, the various specifications of the different sprays and their characteristics are depicted by removing 200 mm from the nozzle exit.

- The round jet occurs without the introduction of pattern air
- The extreme flat spray occurs when the proportion of pattern air (PA) is too high
- The ideal ABC spray occurs when there is a balance between atomising air and pattern air

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
#### Round jet

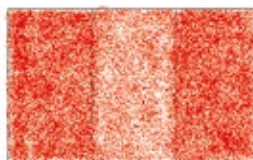Over-humidification in the centre (high flow rate density)  
Finest drops in the area at the edges

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

#### Extreme flat spray

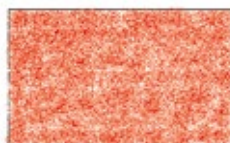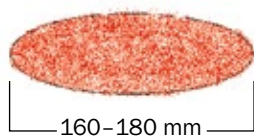Side jets with a high volume current density, overlapping spray patterns lead to the edges being too wet, large droplets, and a fine coverage in the centre

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### PDA measurement technology – measurable success

SCHLICK uses a drop measurement device designed according to the dual PDA principle (Phase-Doppler Anemometry), with a 5-watt (argon-ionic) continuous wave laser.

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#### The right pressure means the right spray

The optimal pressure for atomising and pattern air is controlled by the parameters, quantity of liquid, density and viscosity as well as solid content: You must accommodate the respective application.

Kombinationsmöglichkeiten <i>Combination options</i>	Zerstäubungsluft (AA) in bar (ü) <i>Atomising air (AA) in bar (g)</i>	Formierluft (PA) in bar (ü) <i>Pattern air (PA) in bar (g)</i>
 <b>Feinere Zerstäubung!</b> <b>Fine atomisation!</b>	0.7	0.7 - 1.0 - 1.5
	1.0	1.0 - 1.5 - 2.0
	1.5	1.5 - 2.0 - 2.5 - 3.0
	2.0	2.0 - 2.5 - 3.0 - 3.5
	2.5	2.5 - 3.0 - 3.5

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The values given in the tables refer to the production nozzles S35 an S45.

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The bore diameter of the nozzle should be dependent on the viscosity of the liquid. Here is an example for the production nozzle:

Flüssigkeitsmenge in g/min <i>Quantity of liquid in g/min</i>	Bohrung in mm (Viskosität < 100 mPas) <i>Bore diameter in mm (Viscosity &lt; 100 mPas)</i>	Bohrung in mm (Viskosität > 100 mPas) <i>Bore diameter in mm (Viscosity &gt; 100 mPas)</i>	Min. empfohlener Zerstäubungsluftdruck (AA) in bar (ü) <i>Minimum recommended atomising air (AA) pressure in bar (g)</i>
30 - 60	0.8 - 1.0	1.0 - 1.2	0.7
60 - 120	1.0 - 1.2	1.2 - 1.5	1.0
120 - 150	1.2 - 1.5	1.5 - 1.8	1.5
150 - 180	1.5 - 1.8	1.8 - 2.2	2.0

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The values given in the tables are recommendations for the production nozzles S35 an S45.

Anti-Bearding-Technik<sup>®</sup> speziell für Pharma & Food  
*Anti-Bearding Technology<sup>®</sup> especially for Pharma & Food*

ABC/PCA-Technik (Pharma & Food)  
ABC/PCA Technique (Pharma & Food)

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- Coating
- Befeuchten
- Beschichten
- Beleimen
- Benetzen

## Applications

- Coating
- Humidifying
- Treating
- Gluing
- Moistening

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*The nozzle for production facilities with combined pattern air (PA) and atomising air (AA) regulation provides facilities with the advantage of ABC technology with only one air supply*

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*ABC technology with separate pattern air (PA) and atomising air (AA) regulation, especially for production facilities*

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*Perfect upscaling enables users to proceed from table-top devices to entire laboratory systems, up to and including production itself – all with the aid of the SCHLICK ABC Technique®.*

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*ABC technology with separate pattern air (PA) and atomising air (AA) regulation, especially for laboratory facilities*

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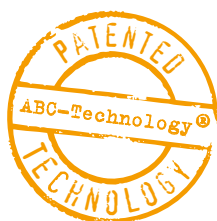
*Downscaling for perfect production processes*

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
*Use perfect nozzle technology right from the smallest batch size! Our latest development, the SCHLICK NANO ABC nozzle, combines all the features of patented SCHLICK ABC technology. Specially developed for atomising the smallest quantities, the NANO nozzle is perfect for conducting initial tests on table-top devices.*

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*Low maintenance requirements ensure considerable cost savings. Thanks to the patented SCHLICK Anti-Bearding Cap (ABC), caking and product build-up are largely eliminated.*

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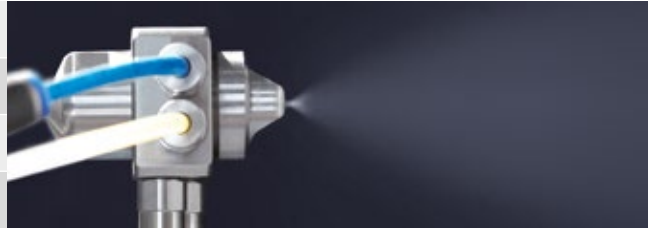*A special stop position attaches the air cap, ensuring results remain uniform and operational errors are eliminated.*

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*GMP-compliant design: extremely easy assembly and disassembly for cleaning and maintenance mean less downtime and more efficient processes.*

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
*Separate connections for pattern air (PA) and atomising air (AA) allow more freedom for adjustment. The atomising air connection can also be used as a retaining tube (M5). Min ID = 2.0 mm. The nozzle is also available with liquid return.*

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*Downscaling for perfect production processes: The SCHLICK NANO ABC nozzle has been specially developed for small drums (8.5", 10.5" and 12"), making the SCHLICK ABC technique now available for very small systems.*

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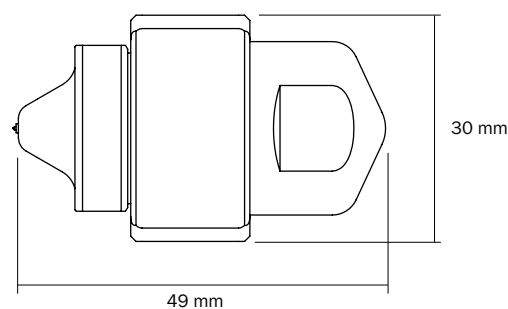Test Center  
Test Center

# Modell 970 Form 7-1 S75 Model 970 Design 7-1 S75

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*Better performance  
for production facilities*

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*GMP design: Very simple installation/de-installation for cleaning or servicing by hand. This leads to shorter downtimes and more efficiency in production.*

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*Cost savings through significant reductions in servicing. Thanks to the SCHLICK patented Anti-Bearding Cap (ABC) the sticking and build-up of deposits is prevented as far as possible.*

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*Standard bore holes for liquid operation: 0.5/0.8/1.0/1.2 mm. Alternative bore hole diameters available on request.*

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*For the first time upscaling from laboratory to production is possible due to the use of ABC-Technology® throughout.*

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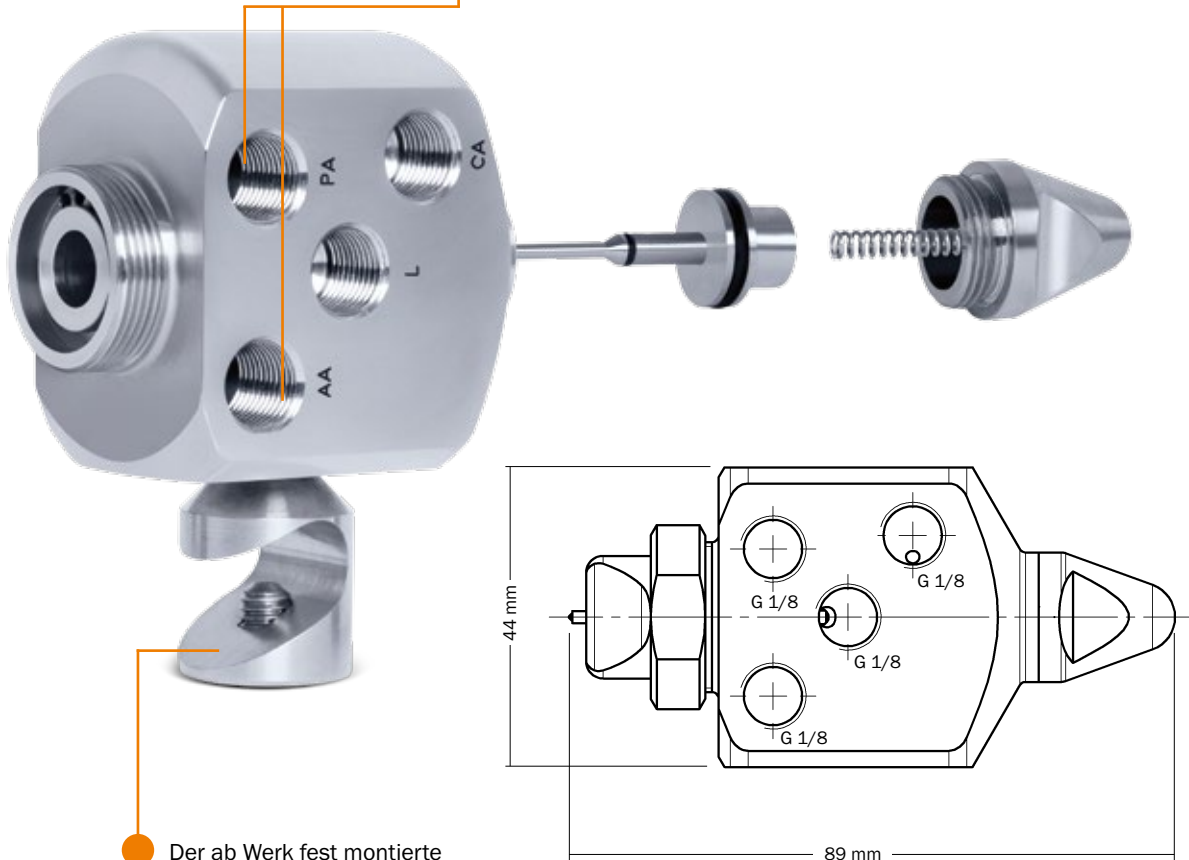
*All attachments are on one side. This leads to a better overview and fewer operator errors.*

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*Separate attachments for pattern air (PA) and atomising air (AA) allow more freedom for calibration. This can help better compensate for turbulence in smaller drums. Available as a model with return system attachment or Design O without needle control (see “custom designs”).*

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*The factory fitted mounting block allows various options for installation.*

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

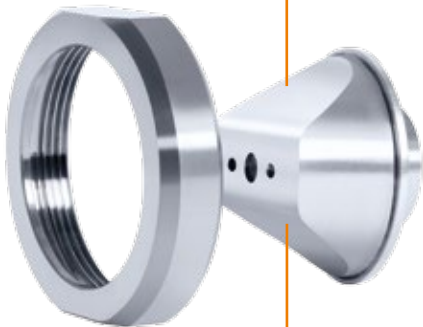
*Standard bore holes for liquid operation: 0.8/1.2/1.5/1.8/2.2 mm. Alternative bore hole diameters available on request.*

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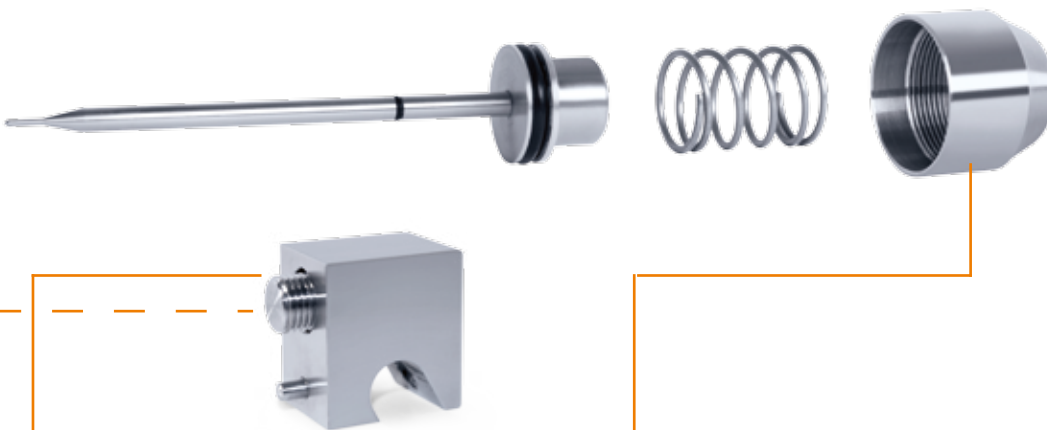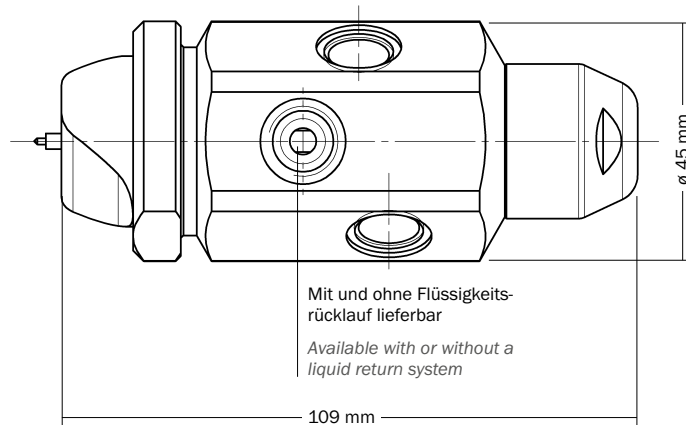*Through separate attachments for the pattern air (PA) and the atomising air (AA) there is more freedom for adjustment. Using the air pressure there are equally as simple and reproducible settings for the droplet size and the spray angle. Available with a liquid return system.*

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*Fixing unit for attachment to a rod available separately. Standard diameter 10 mm – other rod diameters available on request.*

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Better performance  
for production facilities

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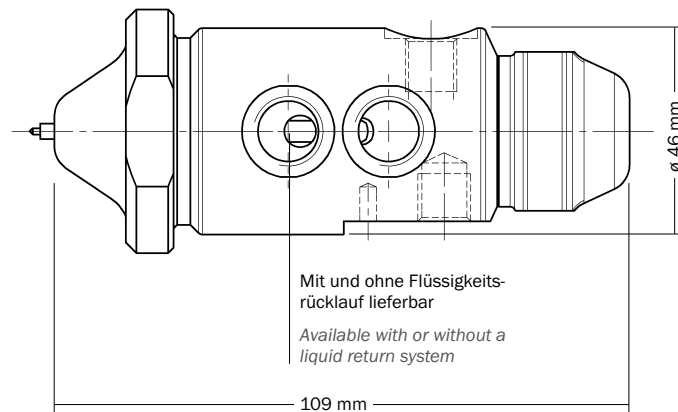
*The spraying angle is set through the various slats in the body of the nozzle.*

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*A single attachment for pattern air (PA) and atomising air (AA). This allows facilities with only one air supply to benefit from the advantage of ABC technology. Available with a liquid return system.*

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*GMP design: Very easy installation/de-installation by hand for cleaning or servicing allows for shorter downtimes and more efficiency in production.*

## SCHLICK Professionell Coating Arm – Der Coating-Spezialist *SCHLICK Professional Coating Arm – The Coating specialist*

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*The SCHLICK PCA is a spray arm especially designed for coating applications. It consists of individual blocks, each with one nozzle. The individual attachment mechanism guarantees use with almost all coating systems.*

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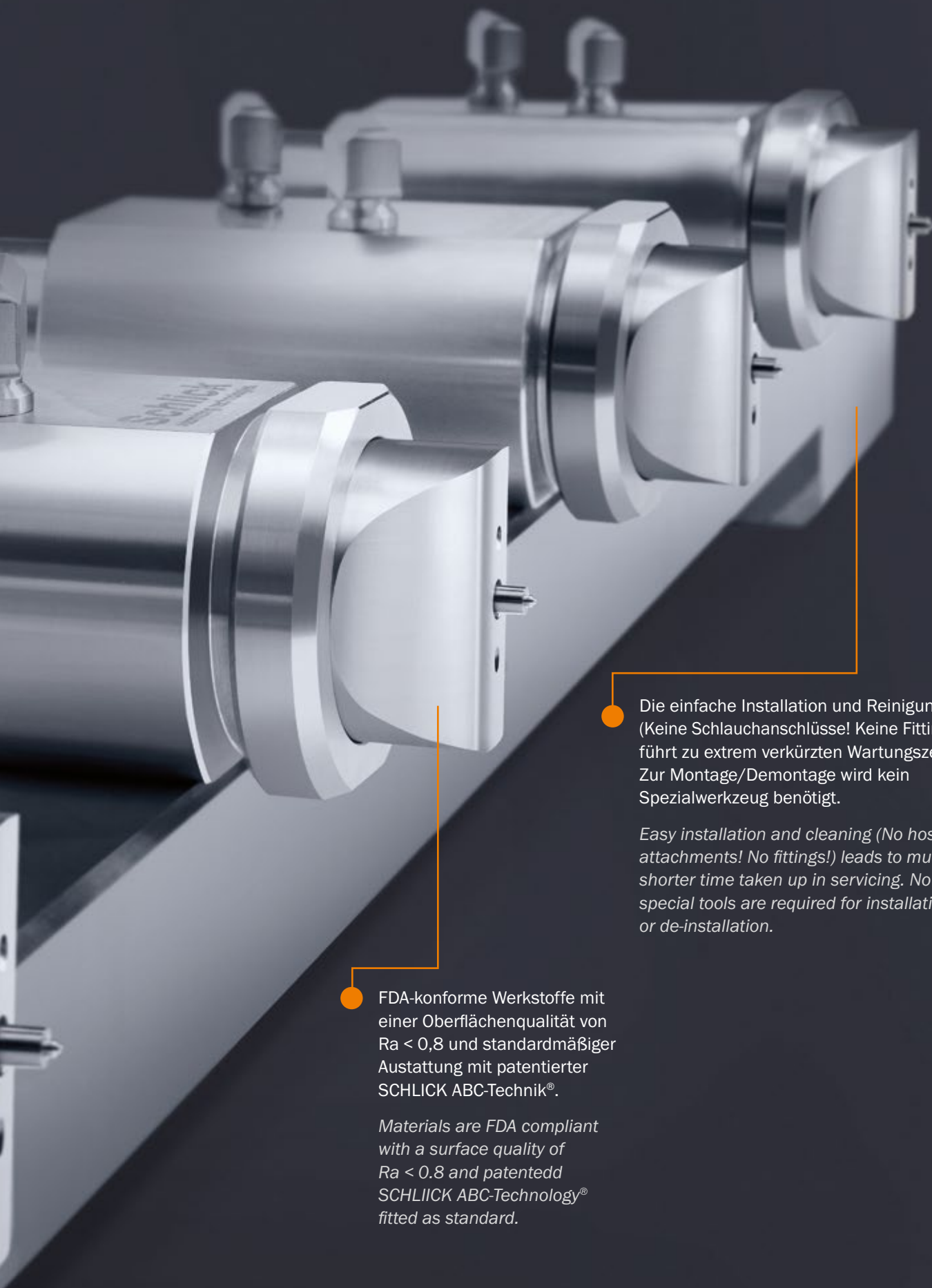
### Materials

*In order to reduce the weight, most of the components are manufactured from titanium 3.7035. The remaining parts are all made from stainless steel 1.4404 (316L). Standard EPDM O-rings (FDA-compliant) are used.*

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*Materials are FDA compliant with a surface quality of  $Ra < 0.8$  and patented SCHLICK ABC-Technology® fitted as standard.*

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*Easy installation and cleaning (No hose attachments! No fittings!) leads to much shorter time taken up in servicing. No special tools are required for installation or de-installation.*

- Produktübersicht  
Product Overview
- Zweistoffdüsen  
Two-Substance  
Nozzles
- ABC/PCA-1 Technik  
ABC/PCA Technique
- Drei-/Vier-Stoffdüsen  
Three-/Four-  
Substance Nozzles
- Multispray  
Multispray
- Einsteckrohre  
Insertion Pipes
- Düsenköpfe  
Nozzle Heads
- Vollkegeldüsen  
Full-Cone Nozzles
- Hohlkegeldüsen  
Hollow-Cone Nozzles
- Hartmetaldüsen  
Carbide Nozzles
- Flachstrahldüsen  
Flat-Spray Nozzles
- Glattdüsen  
Smooth-Jet Nozzles
- Mischdüsen  
Mixing Nozzles
- Dampf-/  
Luftstrahldüsen  
Laval Nozzles
- Reinigungsdüsen  
Cleaning Nozzles
- Zubehör  
Accessories
- Test Center  
Test Center

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*Better performance  
for coating-processes*

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*SCHLICK model 930 Design 7-1 S37: Custom models of the proven S35 model for mounting on the PCA with only two screws. All media are fed through the specially developed PCA connection. There are no further connectors; dirty corners are avoided.*

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### Clever SCHLICK technology

No special tools are required for installation or de-installation – only two screw fittings need to be loosened. The individual nozzles are only fixed with two screws. One individual nozzle in turn consists of only seven individual pieces and O rings.

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*The individual attachment mechanism guarantees use with all coating systems.*

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*The construction system allows the variable use of two to eight nozzles. Advantage: Errors in separation distance between the nozzles arising through installation are eliminated through clearly defined dimensions of the blocks.*

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*Delivery includes individual control air, atomising air, pattern air attachments and liquid attachment with supply and return system.*

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## SCHLICK PCA – The Coating Specialist in Detail

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### Clever technology for coating

The SCHLICK PCA (Professional Coating Arm) is a spraying arm developed **especially for coating applications**. It consists of individual blocks, each with one nozzle.

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The construction system allows the variable use of two to eight nozzles. **Advantage: Through clearly defined dimensions of the blocks, errors in separation distance between the nozzles arising through installation are eliminated.**

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The attachment mechanism for the PCA is individually adapted to the relevant coater.

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### PCA plus ABC: the ideal combination

The PCA nozzles are **fitted with the new SCHLICK anti-bearding technology (ABC)** and a cleaning needle as standard.

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Depending on design the individual nozzles do not require hose attachments or fittings. This completely avoids dirty corners and dramatically reduces cleaning times.

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*Perfection to the last detail*

*Take advantage of our comprehensive expertise – from design to installation – the conception of new products or the optimisation of existing plant. We would be glad to help you improve the success of your operation.*

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*PCA spraying pattern with two-substance nozzles*

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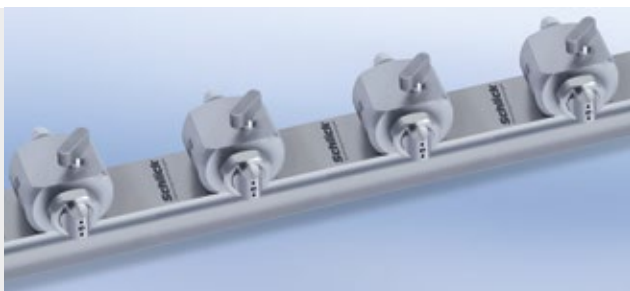
*PCA spraying pattern with pressure nozzles*

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*Due to the substitution of the two-substance nozzles in comparison with the flat spray pressure nozzles, more simple and quick conversion to a **sugar-coating arm** is possible.*

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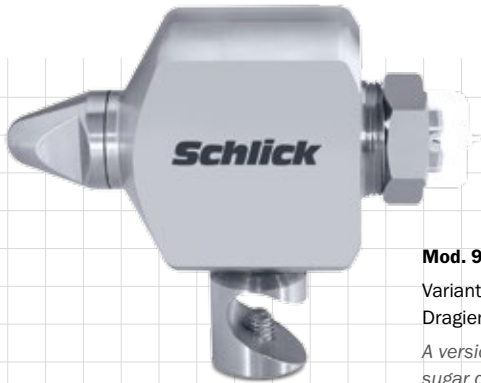
### ***Custom and individual designs, small series***

*As a leading nozzle manufacturer, high quality standardised solutions are business-as-usual for SCHLICK.*

*However, thanks to the breadth of its manufacturing capacities and comprehensive expertise, SCHLICK can also produce custom and individual designs for specific applications.*


*No matter how small the series or short the timeframe.*

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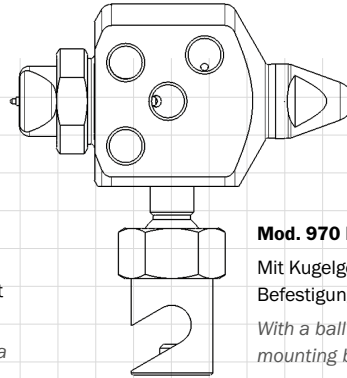
*A version of model 970 S75 with a sugar coating modification unit.*

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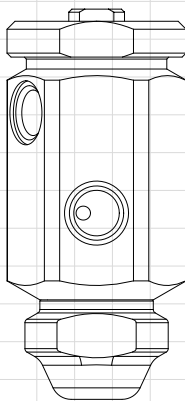
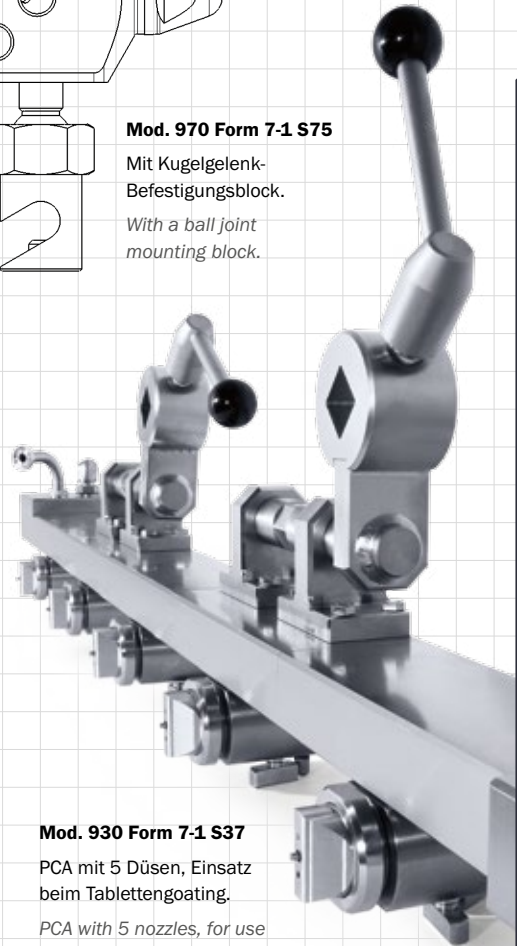
*A custom design of model 930 S75 with a sugar coating modification unit.*

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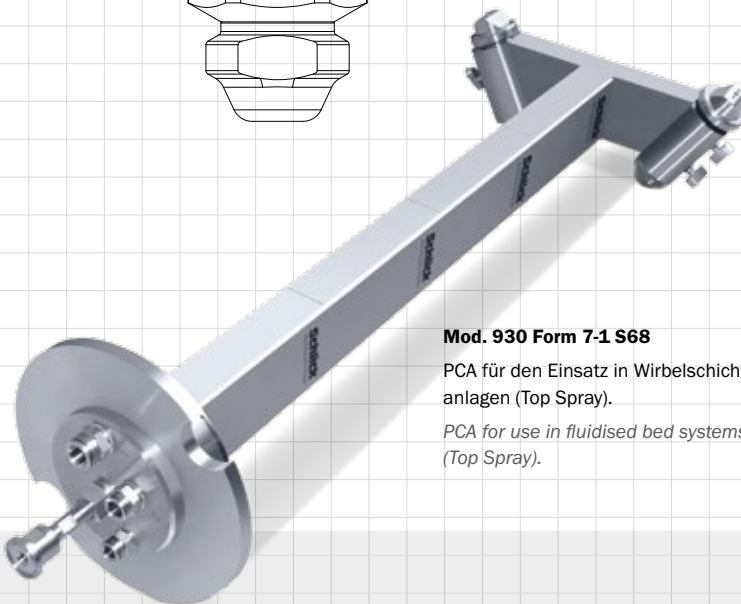
*PCA with 5 nozzles, for use in tablet coating.*

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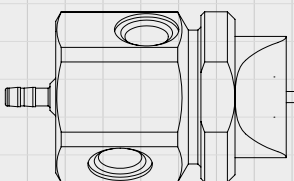
*PCA for use in fluidised bed systems (Top Spray).*

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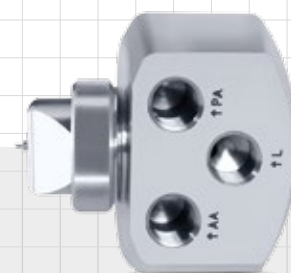*Significantly simplified design with special liquid insert. Absolutely no deviation of the flow direction. For highly sticky and problematic liquids.*

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*Simplified model without needle control. Available with and without mounting block.*

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